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Sciences and Medical Studies
Laboratories Affairs Department
Lab Safety Unit
LS-SciMed@ksu.edu.sa 011 8050052

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What you should do if you need to have Picric Acid in your lab:

1. Make sure that the Picric Acid is kept wet.
2. Do not open a new bottle until needed. Then date the container to show when it was first used to help you in a routine inspection program.
3. Check the hydration water as necessary.
4. Do not use metal spatulas to remove the material.
5. Be sure to clean the bottleneck, cap and threads with a wet cloth before resealing.
6. Get rid of old bottles with metal caps.
7. Do not store large amounts of Picric Acid. Dispose of your Picric Acid every two years.

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When handling Picric Acid:

1. Avoid direct contact - wear appropriate personal protective equipment (Eye Protection, safety glasses, closed lab coat with sleeves fully extended to the wrist lab coat, Gloves)
2. Work in a chemical fume hood when dispensing Picric Acid.
3. Do not break the seal on the container until the material is needed.
4. Visually inspect the container prior to use.
5. Do not use metal spatulas to remove the material from the container.
6. Before closing a container of Picric Acid, always clean the neck of the bottle, cap and thread with a wet cloth before recapping and seal cap with Parafilm.
7. Do not allow Picric Acid to come into contact with metal that is readily oxidized, and do not transfer the material to a container with a metal cap.
8. Never work alone when using Picric Acid.

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Before handling Picric Acid:

1. Read the Material Safety Data for Picric Acid.
2. Identify work practices, engineering controls and protective equipment necessary to do the experiment safely, and include them in written lab procedures.
3. Ensure that engineering controls are operating and protective equipment is available.
4. Go through a dry run to identify potential hazards before using Picric Acid.
5. Be prepared for accidents - know the location of all safety equipment, including eyewash and the first aid kit, emergency contacts and phone numbers, and keep your coworkers informed about your activities with Picric Acid.
6. Do not assume that all lab personnel know the hazards associated with the use of Picric Acid.

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